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Tube or Bottle	Filling Level Requirements	
	Swinging Bucket Rotors	Fixed Angle Rotors
Polyallomer thinwall tubes thickwall tubes Quick-Seal tubes bottles	within 2 to 3 mm of top at least 1/2 full full and heat sealed min to max (see rotor manual) with screw-on cap or cap assembly	full with cap 1/2 full to max capless level or full with cap full and heat sealed 1/2 full to max (see rotor manual) with screw-on cap or cap assembly
Ultra-Clear open-top tubes Quick-Seal tubes	within 2 to 3 mm of top not used	full with cap full and heat sealed
Polycarbonate thickwall tubes bottles	at least 1/2 full at least 1/2 full	1/2 full to max capless level or full with cap min to max (see rotor manual) with screw-on cap or cap assembly
Stainless Steel tubes	any level	any level with cap or cap assembly
Polypropylene tubes and bottles	at least 1/2 full	1/2 full to max capless level or full with cap or cap assembly
Polyethylene tubes	at least 1/2 full	1/2 full to max capless level or full with cap
Cellulose Propionate tubes and bottles	at least 1/2 full	1/2 full to max capless level

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