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Tube or Bottle	Filling Level Requirements	
	Swinging Bucket Rotors	Fixed Angle Rotors
<b>Polyallomer</b> thinwall tubes thickwall tubes  Quick-Seal tubes bottles	within 2 to 3 mm of top at least 1/2 full  full and heat sealed min to max (see rotor manual) with screw-on cap or cap assembly	full with cap 1/2 full to max capless level or full with cap full and heat sealed 1/2 full to max (see rotor manual) with screw-on cap or cap assembly
<b>Ultra-Clear</b> open-top tubes Quick-Seal tubes	within 2 to 3 mm of top not used	full with cap full and heat sealed
<b>Polycarbonate</b> thickwall tubes  bottles	at least 1/2 full  at least 1/2 full	1/2 full to max capless level or full with cap min to max (see rotor manual) with screw-on cap or cap assembly
<b>Stainless Steel</b> tubes	any level	any level with cap or cap assembly
<b>Polypropylene</b> tubes and bottles	at least 1/2 full	1/2 full to max capless level or full with cap or cap assembly
<b>Polyethylene</b> tubes	at least 1/2 full	1/2 full to max capless level or full with cap
<b>Cellulose Propionate</b> tubes and bottles	at least 1/2 full	1/2 full to max capless level

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